

MassFinder 3 Keyboard Shortcuts

You can control MassFinder with the mouse. Only administrators or advanced users have to use keyboard commands. The need to type specific keyboard commands protects regular users from accidentally perform serious modification. Anyway, make sure to deactivate *allow library modification* at all times (except when actually administrating the library).

+	add library entry
* (asterisk)	view/modify library entry
BACKSPACE	delete library entry
SPACE	adjust current GC/MS scan to peak (jump to nearest maximum)
n	edit name of current GC/MS scan (selection box)
N	edit name of current GC/MS scan (simple method)
E	edit library spectrum
!	jump to GC/MS scan by number
' (APOSTROPH)	jump to library entry by internal number
=	jump to library entry by unique key
#	jump to library entry by ordered list index
<	jump to previous marked GC/MS peak
>	jump to next marked GC/MS peak
e	toggle on/off enhanced search algorithm
l (lower-case L)	toggle on/off library-based comparison
r	toggle on/off two-dimensional search algorithm (use of retention indices)
s	toggle on/off alternative subtraction layout

(save GC/MS (with labelled peaks and retention indices)
)	save library
b	show baseline on/off
q	show quality hint on/off
X	remove all peak labels of GC/MS
M	mark library entry as "manual"
A	mark library entry as "administrator"